

AmeriDx® PLANT DNA QUICK & EASY PURIFICATION KIT



Easy to use, Two Steps and
Easy to Scale up

Quick & Easy DNA Purification Kit


Unlock the future of plant DNA extraction with our Plant DNA Quick & Easy Purification Kit! This groundbreaking kit offers a seamless two-step process—lysis and purification—that ensures the highest quality DNA ready for downstream applications like PCR and qPCR. No equipment required means you can streamline your workflow with ease, scaling up to high-throughput with consistent, reliable results. Say goodbye to complex protocols and hello to efficient, superior DNA purification. Get yours today and revolutionize your research!

Assays

1. Cut to several pieces of the plant leaves/root/stem, and mix with Lysis buffer, incubate at -85°C for 20 minutes
2. Take the solution to another Eppendorf tube, mix with equal volume of Extraction Buffer
3. Spin at 12,000 rpm for 5 minutes, save the supernatant for downstream assays (PCR, qPCR etc).

Sensitivity and Specificity

Work with 5 mm plant tissues, retrieve >100 uL DNA solution.

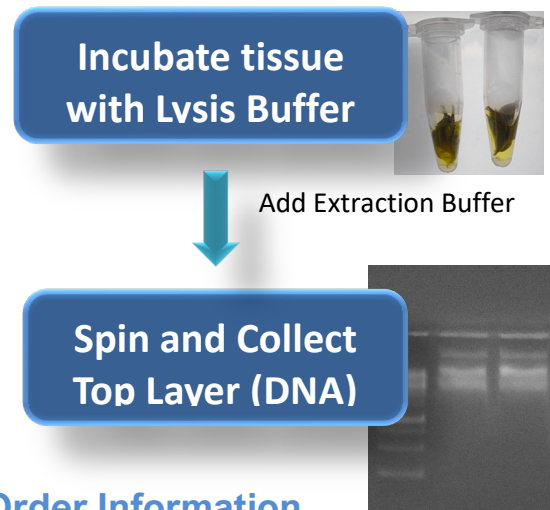

Order Information

Product Name	Catalog No
Plant DNA Quick&Easy Purification Kit	Rxxxx
Microbe DNA Quick&Easy Purification Kit	Rxxxx



ivena Ltd. Pure Offices,
to Close, Warwick CV34 6WE UK



AmeriDx Diagnostics, Inc

9235 Activity Road Suite 101, San Diego, CA 92126 USA

Tel: +1 (858) 282-8567 | Info@ameridx.com | <https://www.ameridx.com>